

Fibronectin IHC-Fr

Fix slides in ice-cold acetone at -20°C for 5 min

Rehydrate 2X 5 minutes in 1X PBS at RT

Block for 1 hour minutes with 10% donkey serum in 1X PBS in humid chamber at RT

Incubate for 2 hours with Rabbit Anti Fibronectin (1:50 dilution) in 10% donkey serum in 1X PBS in humid chamber at RT

Wash 2X 5 minutes in 1X PBS

Incubate for 2 hours with Rhodamine Red X Donkey Anti Rabbit IgG (1:50 dilution) in 10% donkey serum in 1X PBS in humid chamber at RT

Wash 2X 5 minutes in 1X PBS in a covered container

Rinse briefly in d H2O and coverslip with 1-2 drops of selected mounting medium